




DONNING GLOVES

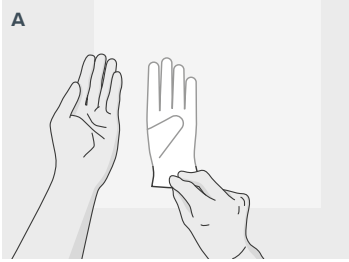


Grasp the interior of one glove and carefully pull over the opposite hand

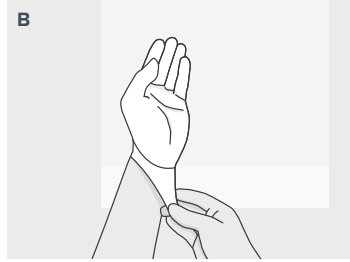


Gently ease the glove up over the wrist to cover gown sleeve

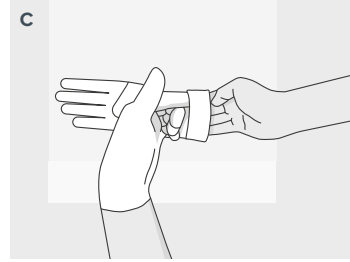


Lift the second glove under the folded cuff using a gloved finger

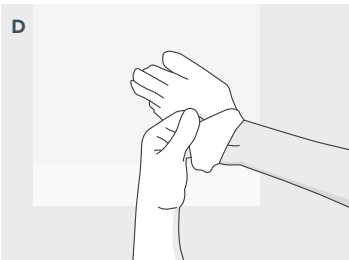

Gently ease your second hand into the other glove

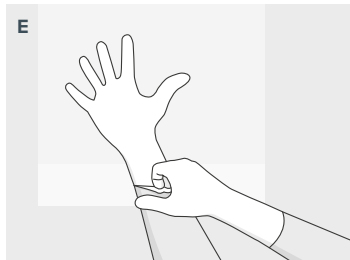


Using a gloved finger, unfold the cuff over your wrist and gown sleeve

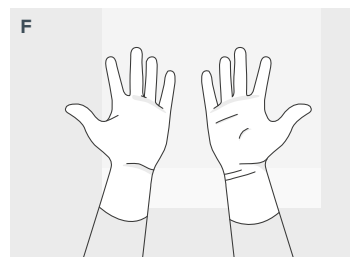


Make sure your gloved fingertips do not touch bare forearms or wrists


DOFFING GLOVES

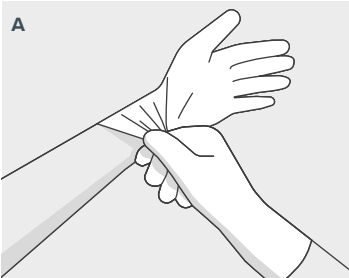


Grasp the exterior of one glove with your other gloved hand.

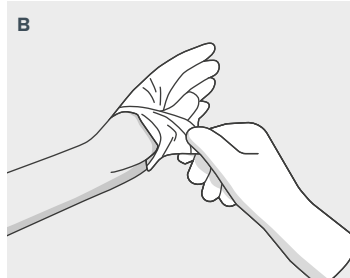


Carefully pull the glove off your hand, turning it inside-out.

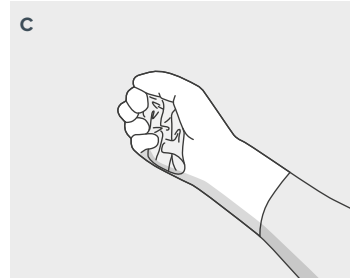

Ball the glove up and hold in your other gloved hand.

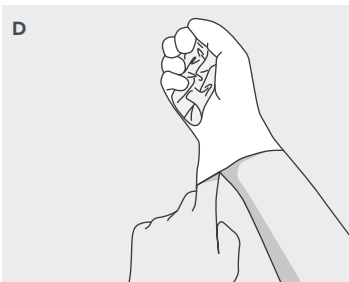

Slide your ungloved finger into the opening of the other glove.

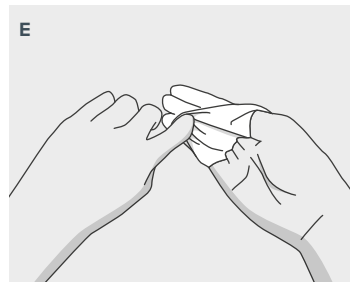

Carefully pull the glove off your hand, turning it inside out again.



Discard appropriately.

Distributed by Fisher Scientific. Contact us today:

Austria: fishersci.at Belgium: fishersci.be Denmark: fishersci.dk
Germany: fishersci.de Ireland: fishersci.ie Italy: fishersci.it
Finland: fishersci.fi France: fishersci.fr Netherlands: fishersci.nl
Norway: fishersci.no Portugal: fishersci.pt Spain: fishersci.es
Sweden: fishersci.se Switzerland: fishersci.ch UK: fishersci.co.uk

© 2022 Thermo Fisher Scientific Inc. All rights reserved.
Trademarks used are owned as indicated at fishersci.com/trademarks.