

KIMTECH™

Kimtech™ G3 Sterile Sterling™ Nitrile Gloves

Donning Technique

Before starting the donning procedure, wash hands thoroughly and dry.

1

Peel open sterile pouch and unfold glove wallet (DO NOT touch the exterior surface of gloves). Pinch the sides of wallet to open.



2

Apply first glove to hand by sliding palm up into glove (thumb facing outward). Bend thumb toward center of palm and slide into glove while pulling up on the cuff. Leave the cuff rolled up.



3

Apply second glove to hand by sliding the four gloved fingers into cuff of the second glove. Slide ungloved palm (thumb facing outward) into glove. Bend thumb toward center of palm and slide into glove while pulling up with fingers of gloved hand.



4

Complete donning the gloves by pulling up the cuff of the first glove with the fingers of the second hand.



5

If double donning is desired, repeat steps 1-4 with a second set of gloves - this time using a half inch size larger glove if necessary.



Visit our website at www.kimtech.eu
Technical help: kimtech@kcc.com

©/™ Trademarks of Kimberly-Clark Worldwide, Inc. ©2020 KCWW Publication code ID 5372.01 EN 06.20

Distributed by Fisher Scientific. Contact us today:

Austria: fishersci.at Belgium: fishersci.be Denmark: fishersci.dk
Germany: fishersci.de Ireland: fishersci.ie Italy: fishersci.it
Finland: fishersci.fi France: fishersci.fr Netherlands: fishersci.nl
Norway: fishersci.no Portugal: fishersci.pt Spain: fishersci.es
Sweden: fishersci.se Switzerland: fishersci.ch UK: fishersci.co.uk

© 2022 Thermo Fisher Scientific Inc. All rights reserved.
Trademarks used are owned as indicated at fishersci.com/trademarks.

 **fisher scientific**
part of Thermo Fisher Scientific